



KERN & Sohn GmbH

Ziegelei 1
D-72336 Balingen
E-mail:
info@kern-sohn.com

Tel.: +49-[0]7433-9933-0
Fax: +49-[0]7433-9933-149
Internet:
www.kern-sohn.com

Operating Manual Precision balance

KERN PFB

Version 1.7
01/2012
GB



PFB-BA-e-1217



KERN PFB

Version 1.7 01/2012

Operating Manual

Precision balance

Table of Contents

1	Technical Data	4
2	Declaration of conformity	6
3	Basic Information (General)	7
3.1	Proper use	7
3.2	Improper Use	7
3.3	Warranty	7
3.4	Monitoring of Test Resources.....	8
4	Basic Safety Precautions	8
4.1	Pay attention to the instructions in the Operation Manual.....	8
4.2	Personnel training.....	8
5	Transportation & Storage	8
5.1	Testing upon acceptance	8
5.2	Packaging / return transport	8
6	Unpacking, Setup and Commissioning	9
6.1	Installation Site, Location of Use	9
6.2	Unpacking.....	9
6.2.1	Placing/ Remove the transportation lock.....	10
6.2.2	Scope of delivery	10
6.3	Mains connection.....	10
6.4	Connection of peripheral devices	10
6.5	Initial Commissioning.....	11
6.6	Adjustment.....	11
6.7	Adjusting.....	11
6.8	Linearization	13
6.8.1	Table Linearization points.....	15
7	Controls	16
7.1	Keyboard overview	16
7.2	Overview of display	17
8	Operation	18
8.1	Simple weighing	18
8.2	Taring.....	19
8.3	Parts counting.....	20
8.4	Adding.....	21
8.5	Percent weighings	23

9	The menu	24
9.1	Menu map.....	24
9.2	Call up menu	24
9.3	Navigation in the menu.....	24
9.4	Menu overview	24
9.5	Menu settings	26
9.5.1	Switch over weighing units (Unit)	26
9.5.2	Display background illumination.....	27
10	Data output RS232	28
10.1	Technical data	28
10.2	Pin allocation of balance output bushing:.....	28
10.3	Printer operation - printout examples	29
10.4	Continuous data output	29
10.5	Remote control instructions	30
11	Error messages.....	30
12	Service, maintenance, disposal.....	31
12.1	Cleaning.....	31
12.2	Service, maintenance	31
12.3	Disposal	31
13	Instant help.....	32

1 Technical Data

KERN	PFB 120-3	PFB 200-3	PFB 300-3	PFB 1200-2
Readability (d)	0,001 g	0,001 g	0,001 g	0,01 g
Weighing range (max)	120 g	200 g	300 g	1200 g
Reproducibility	0,001 g	0,001 g	0,001g	0,01 g
Linearity	±0,003 g	±0,003 g	±0,003 g	±0,03 g
Smallest part weight for piece counting	2 mg	2 mg	2 mg	20 mg
Reference quantities at piece counting	10, 20, 50, 100, 200			
Weighing Units	g, ct, lb, oz, d, ozt, dwt, mo, tl h, tl c, tl t, t, bt, n			
Recommended adjustment weight, not added (class)	100 g (F1)	200 g (F1)	300 g (F1)	1000 g (F1)
Warm-up time	2 hours			
Stabilization time (typical)	3 sec.			
Operating temperature	+ 5° C + 35° C			
Humidity of air	max. 80 % (not condensing)			
Housing (B x D x H) mm	200 x 257 x 152 (with draft shield) 200 x 257 x 87 (without draft shield)			
Draft shield (B x D x H) mm, plastic	158 x 143 x 64 (inside) 167 x 154 x 80 (outside)			
Weighing plate, stainless steel (mm)	Ø 80	Ø 80	Ø 80	Ø 120
Weight kg (net)	2 kg			
Underfloor weighing	optional			
Voltage	220-240 V / 50 Hz			

KERN	PFB 2000-2	PFB 3000-2	PFB 6000-1	PFB 6K0.05
Readability (d)	0,01 g	0,01 g	0,1 g	0,05 g
Weighing range (max)	2000 g	3000 g	6000 g	6000 g
Reproducibility	0,01 g	0,01 g	0,1 g	0,05 g
Linearity	±0,03 g	±0,03 g	± 0,3 g	± 0,15 g
Smallest part weight for piece counting	20 mg	20 mg	200 mg	200 mg
Reference quantities at piece counting	10, 20, 50, 100, 200			
Weighing Units	g, ct, lb, oz, d, ozt, dwt, mo, tl h, tl c, tl t, t, bt, n			
Recommended adjustment weight, not added (class)	2000 g (F1)	3000 g (F1)	5000 g (F1)	5000 g (F1)
Warm-up time	2 hours			
Stabilization time (typical)	3 sec.			
Operating temperature	+ 5° C + 35° C			
Humidity of air	max. 80 % (not condensing)			
Housing (B x D x H) mm	200 x 257 x 152 (with draft shield) 200 x 257 x 87 (without draft shield)		-	-
Draft shield (B x D x H) mm, plastic	158 x 143 x 64 (innen) 167 x 154 x 80 (außen)		-	-
Weighing plate, stainless steel (mm)	Ø 120	Ø 120	155 x 145	155 x 145
Weight kg (net)	2 kg			
Underfloor weighing	optional			
Voltage	220-240 V / 50 Hz			

2 Declaration of conformity



KERN & Sohn GmbH
 D-72322 Balingen-Frommern
 Postbox 4052
 email: info@kern-sohn.de

Phone: 0049-[0]7433- 9933-0
 Fax: 0049-[0]7433-9933-149
 Internet: www.kern-sohn.de

Declaration of conformity

EC-Konformitätserklärung
EC- Déclaration de conformité
EC-Dichiarazione di conformità
EC- Declaração de conformidade
EC-Deklaracja zgodności

EC-Declaration of -Conformity
EC-Declaración de Conformidad
EC-Conformiteitverklaring
EC- Prohlášení o shode
ЕС-Заявление о соответствии

D	Konformitäts- erklärung	Wir erklären hiermit, dass das Produkt, auf das sich diese Erklärung bezieht, mit den nachstehenden Normen übereinstimmt.
GB	Declaration of conformity	We hereby declare that the product to which this declaration refers conforms to the following standards.
CZ	Prohlášení o shode	Tímto prohlašujeme, že výrobek, kterého se toto prohlášení týká, je v souladu s níže uvedenými normami.
E	Declaración de conformidad	Manifetamos en la presente que el producto al que se refiere esta declaración está de acuerdo con las normas siguientes
F	Déclaration de conformité	Nous déclarons avec cela responsabilité que le produit, auquel se rapporte la présente déclaration, est conforme aux normes citées ci-après.
I	Dichiarazione di conformità	Dichiariamo con ciò che il prodotto al quale la presente dichiarazione si riferisce è conforme alle norme di seguito citate.
NL	Conformiteit- verklaring	Wij verklaren hiermede dat het product, waarop deze verklaring betrekking heeft, met de hierna vermelde normen overeenstemt.
P	Declaração de conformidade	Declaramos por meio da presente que o produto no qual se refere esta declaração, corresponde às normas seguintes.
PL	Deklaracja zgodności	Niniejszym oświadczamy, że produkt, którego niniejsze oświadczenie dotyczy, jest zgodny z poniższymi normami.
RUS	Заявление о соответствии	Мы заявляем, что продукт, к которому относится данная декларация, соответствует перечисленным ниже нормам.

Electronic Balance: KERN PFB

Mark applied	EU Directive	Standards
CE	2004/108/EC EMC	EN 61000-6-1 :2007 EN 61000-6-3 :2007 EN 61000-3-3 : 1995+A1 :2001+A2 :2005 EN 61000-3-2 :2006

Date: 18.09.2008

Signature:

KERN & Sohn GmbH
Management

KERN & Sohn GmbH, Ziegelei 1, D-72336 Balingen, Tel. +49-[0]7433/9933-0, Fax +49-[0]7433/9933-149

English

3 Basic Information (General)

3.1 Proper use

The balance you purchased is intended to determine the weighing value of material to be weighed. It is intended to be used as a “non-automatic” balance, i.e. the material to be weighed is manually and carefully placed in the centre of the weighing plate. As soon as a stable weighing value is reached the weighing value can be read.

3.2 Improper Use

Do not use balance for dynamic weighing. In the event that small quantities are removed or added to the material to be weighed, incorrect weighing results can be displayed due to the “stability compensation” in the balance. (Example: Slowly draining fluids from a container on the balance.)

Do not leave permanent load on the weighing plate. This may damage the measuring system.

Impacts and overloading exceeding the stated maximum load (max) of the balance, minus a possibly existing tare load, must be strictly avoided. Balance may be damaged by this.

Never operate balance in explosive environment. The serial version is not explosion protected.

The structure of the balance may not be modified. This may lead to incorrect weighing results, safety-related faults and destruction of the balance.

The balance may only be used according to the described conditions. Other areas of use must be released by KERN in writing.

3.3 Warranty

Warranty claims shall be voided in case

- Our conditions in the operation manual are ignored
- The appliance is used outside the described uses
- The appliance is modified or opened
- Mechanical damage and damage caused by media, liquids
Natural wear and tear
- The appliance is improperly set up or incorrectly electrically connected
- The measuring system is overloaded

3.4 Monitoring of Test Resources

In the framework of quality assurance the measuring-related properties of the balance and, if applicable, the testing weight, must be checked regularly. The responsible user must define a suitable interval as well as type and scope of this test. Information is available on KERN's home page (www.kern-sohn.com) with regard to the monitoring of balance test substances and the test weights required for this. In KERN's accredited DKD calibration laboratory test weights and balances may be calibrated (return to the national standard) fast and at moderate cost.

4 Basic Safety Precautions

4.1 Pay attention to the instructions in the Operation Manual

Carefully read this operation manual before setup and commissioning, even if you are already familiar with KERN balances.

4.2 Personnel training

The appliance may only be operated and maintained by trained personnel.

5 Transportation & Storage

5.1 Testing upon acceptance

When receiving the appliance, please check packaging immediately, and the appliance itself when unpacking for possible visible damage.

5.2 Packaging / return transport



- ⇒ Keep all parts of the original packaging for a possibly required return.
- ⇒ Only use original packaging for returning.
- ⇒ Prior to dispatch disconnect all cables and remove loose/mobile parts.
- ⇒ Reattach possibly supplied transport securing devices.
- ⇒ Secure all parts such as the glass wind screen, the weighing platform, power unit etc. against shifting and damage.

6 Unpacking, Setup and Commissioning

6.1 Installation Site, Location of Use

The balances are designed in a way that reliable weighing results are achieved in common conditions of use.

You will work accurately and fast, if you select the right location for your balance.

Therefore, observe the following for the installation site:

- Place the balance on a firm, level surface;
- Avoid extreme heat as well as temperature fluctuation caused by installing next to a radiator or in the direct sunlight;
- Protect the balance against direct draughts due to open windows and doors;
- Avoid jarring during weighing;
- Protect the balance against high humidity, vapors and dust;
- Do not expose the device to extreme dampness for longer periods of time. Non-permitted condensation (condensation of air humidity on the appliance) may occur if a cold appliance is taken to a considerably warmer environment. In this case, acclimatize the disconnected appliance for ca. 2 hours at room temperature.
- Avoid static charge of goods to be weighed or weighing container.

Major display deviations (incorrect weighing results) may be experienced should electromagnetic fields (e.g. due to mobile phones or radio equipment), static electricity accumulations or instable power supply occur. Change location or remove source of interference.

6.2 Unpacking

Carefully remove the balance from the packaging and setup balance at the intended workstation.

6.2.1 Placing/ Remove the transportation lock

The balance must be installed in a way that the weighing plate is exactly in horizontal position.

Remove the transportation lock [1] (PFB 120-3, PFB 200-3, PFB 300-3):



6.2.2 Scope of delivery

Serial accessories:

- Balance
- Weighing plate
- Mains power supply
- Operating Manual
- Draft shield (only models PFB 120-3, PFB 200-3, PFB 300-3, PFB 1200-2, PFB 2000-2, PFB 3000-2)

6.3 Mains connection

Power is supplied via the external mains adapter. The stated voltage value must be the same as the local voltage.

Only use original KERN mains adapters. Using other makes requires consent by KERN.

6.4 Connection of peripheral devices

Before connecting or disconnecting of additional devices (printer, PC) to the data interface, always disconnect the balance from the power supply.

With your balance, only use accessories and peripheral devices by KERN, as they are ideally tuned to your balance.

6.5 Initial Commissioning

In order to obtain exact results with the electronic balances, your balance must have reached the operating temperature (see warming up time chap. 1). During this warming up time the balance must be connected to the power supply (mains, accumulator or battery).

The accuracy of the balance depends on the local acceleration of gravity. Strictly observe hints in chapter Adjustment.

6.6 Adjustment

As the acceleration value due to gravity is not the same at every location on earth, each balance must be coordinated - in compliance with the underlying physical weighing principle - to the existing acceleration due to gravity at its place of location (only if the balance has not already been adjusted to the location in the factory). This adjustment process must be carried out for the first commissioning, after each change of location as well as in case of fluctuating environment temperature. To receive accurate measuring values it is also recommended to adjust the balance periodically in weighing operation.

6.7 Adjusting

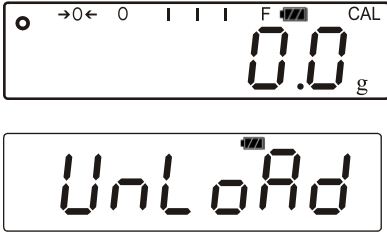


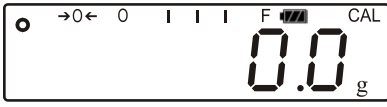
The adjustment should be made with the recommended adjustment weight (see chap. 1 "Technical data").

Procedure when adjusting:

Observe stable environmental conditions.

A warming up time (see chapter 1; Technical Data) is required for stabilization.

Ensure that there are no objects on the weighing plate.

Balance display	Operation
 <p>The display shows a zeroed scale (0.0 g) with indicators for tare (→0←), units (0), and calibration (CAL). Below, the text 'UnLoAd' is displayed.</p>	<p>⇒ In weighing mode press  approx. 3 sec. until „UnLoAd“ appears.</p>
 <p>The display shows the text 'LoAd'.</p>	<p>When display „LoAd“ appears, put the necessary adjustment weight (see chap.1; Technical Data) on the center of the weighing plate.</p> <p>Adjustment will be carried out automatically after dead stop control.</p>
	<p>⇒ Wait for stability display</p>
	<p>„Pass“ appears, the adjustment process has been finished successfully.</p> <p>⇒ Take away adjustment weight</p>
 <p>The display returns to the zeroed scale (0.0 g) with the same indicators as in the first row.</p>	<p>Wait until the balance is again in the weighing mode.</p>

6.8 Linearization

Linearity shows the greatest deviation of a weight display on the scale to the value of the respective test weight according to plus and minus over the entire weighing range.

If linearity deviation is discovered during a testing instrument control, you can improve this by means of linearization.

- In balances with a resolution of > 15 000 dividing steps carrying out a linearization is recommended.
- Carrying out linearization is restricted to specialist staff possessing well acquainted with the workings of weighing scales.
- The test weights to be used must be adapted to the weighing scale's specifications; see chapter "testing instruments control".
- Observe stable environmental conditions. Stabilisation requires a certain warm-up time.
- After successful linearization you will have to carry out calibration; see chapter "testing instruments control".

Balance display	Operation
	⇒ Start balance by pressing Whilst the balance carries out a self test, press until „F1 Unt“ appears
	⇒ Press repeatedly until "tECH" appears
	⇒ Press , „Pin“ appears
	⇒ Press , and successively, „P1 Lin“ appears
	⇒ Press anew, „Pin“ appears again
 	⇒ Press successively , and , „LoAd 0“ appears; after the apparition of the stability display and the signal sound, „LoAd 1“ will be displayed

	<ul style="list-style-type: none"> ⇒ Put on weight 1 (see table chap. 6.8.1) ⇒ Wait for the stability display and the signal sound ⇒ „LoAd 2“ appears ⇒ Remove weight 1 and put on the weight 2
	<ul style="list-style-type: none"> ⇒ After the stability display and the signal sound „LoAd 3“ appears; remove weight 2 and put on weight 3
	<ul style="list-style-type: none"> ⇒ After the stability display and the signal sound „LoAd 4“ appears; remove weight 3 and put on weight 4
	<ul style="list-style-type: none"> ⇒ After having placed weight 4, „LoAd 0“ appears again ⇒ Remove weight 4, „LoAd 4“ appears anew ⇒ Place weight 4 again
	<ul style="list-style-type: none"> ⇒ Wait for the stability display and the signal sound, „LoAd 3“ will be displayed ⇒ Remove weight 4 and place weight 3
	<ul style="list-style-type: none"> ⇒ Wait for the stability display and the signal sound, „LoAd 2“ will be displayed ⇒ Remove weight 3 and place weight 2
	<ul style="list-style-type: none"> ⇒ Wait for the stability display and the signal sound, „LoAd 1“ will be displayed ⇒ Remove weight 2 and place weight 1
	<ul style="list-style-type: none"> ⇒ Wait for the stability display and the signal sound ⇒ „LoAd 0“ will be displayed
	<ul style="list-style-type: none"> ⇒ Remove weight 1
	<ul style="list-style-type: none"> ⇒ After the stability display and the signal sound the balance carries out a selftest and changes into the weighing mode. Now the linearization is concluded successfully.

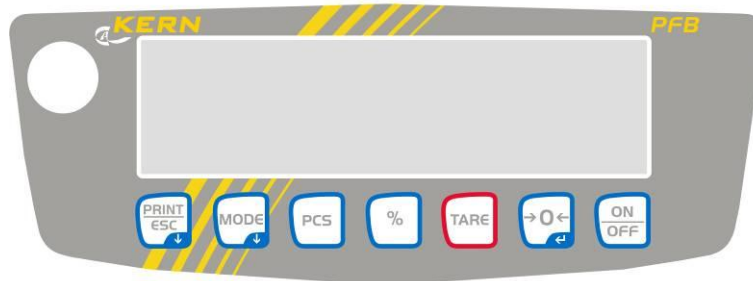
6.8.1 Table Linearization points








Max	Weight 1	Weight 2	Weight 3	Weight 4
120 g	30 g	60 g	90 g	120 g
200 g	50 g	100 g	150 g	200 g
300 g	50 g	100 g	200 g	300 g
1200 g	300 g	600 g	900 g	1200 g
2000 g	500 g	1000 g	1500 g	2000 g
3000 g	0,5 kg	1 kg	2 kg	3 kg
6000 g	1 kg	2 kg	4 kg	6 kg

7 Controls

7.1 Keyboard overview

Buttons specification:



Key	Designation	Pressed once and released
	[ON/OFF]	⇒ Turn on/off
	[ZERO]	<ul style="list-style-type: none"> ⇒ Set weight display at zero ⇒ Activate the current element in the menu ⇒ Select in Weighing units (menu F UNIT): change to the next unit
	[TARE]	⇒ Taring
	[%]	<ul style="list-style-type: none"> ⇒ Call-up %-function ⇒ In the percent mode back to weighing mode ⇒ Only % function
	[PCS]	<ul style="list-style-type: none"> ⇒ Call-up piece counting mode ⇒ In the piece count mode back to weighing mode
	[MODE]	<ul style="list-style-type: none"> ⇒ Weighing units switch-over; ⇒ At weighing unit switch-over (menu „F UNIT“): Set weighing unit on/off ⇒ Counting up in piece counting mode ⇒ Select menu and pass through menu items from top to bottom ⇒ Changing readability
	[PRINT/ESC]	<ul style="list-style-type: none"> ⇒ Print out weighing result ⇒ Exit menu (jumping back to weighing mode)

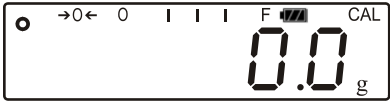

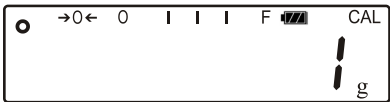
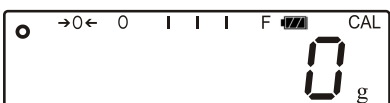

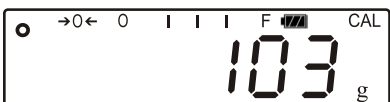


7.2 Overview of display



Display	Description
g	Gram
kg	Kilogram
→0←	Zeroing display
→T←	Taring
o	Stability display
Pcs	Display for piece counting
%	Percent weighing display
◀	Display for tolerance weighing
(mom)	Momme
CAL	Display for adjustment. Signals the adjustment function
0 F	Bar graph
Weight unit display	g (Gramm) (ct) Carat (oz) Ounce (ozt) Fine ounce (d) Dram (dwt) Penny weight (tl.h) Tael (Taiwan) (tl.c) Tael (Chin.) (t.lt) Troy Tael (t) Tola

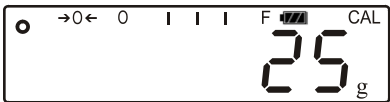
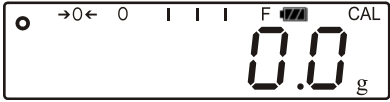

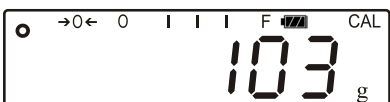
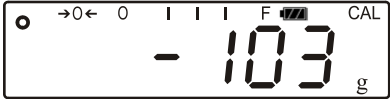
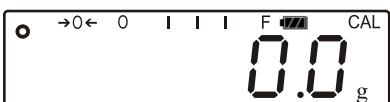

8 Operation

8.1 Simple weighing

Balance display	Operation
 <p>The display shows 0.0 g. Above the display are indicators: a small circle, a zero with left and right arrows, four vertical bars, a battery level indicator (F), and CAL.</p>	<p>⇒ Start balance by pressing  .</p> <p>The balance will carry out a self-test.</p> <p>Wait for „0.0“ display</p>
  <p>The display shows 1 g, then 0 g. The indicators above the display are the same as in the previous row.</p>	<p>⇒ Should the balance not display exactly „0.0“ despite empty scale pan, press the  button.</p> <p>The balance starts with resetting to „0“</p>
 <p>The display shows 103 g. The indicators above the display are the same as in the previous rows.</p>	<p>⇒ Place goods to be weighed on balance</p> <p>Wait until the stability display appears [] .</p> <p>Read weighing result.</p>
	<p>⇒ To switch off the balance press shortly  .</p>

8.2 Taring








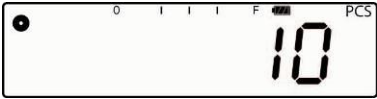




The dead weight of any weighing container may be tared away by pressing a button, so that the following weighings show the net weight of the goods to be weighed.

Balance display	Operation
 <p>The display shows a pictogram of a weighing container, a zero with a left arrow, three vertical bars, a full battery icon, and 'CAL'. The main display shows '25' with a 'g' below it.</p>	<p>⇒ Deposit weighing receptacles. The weight of the container is displayed.</p>
 <p>The display shows a pictogram of a weighing container, a zero with a left arrow, three vertical bars, a full battery icon, and 'CAL'. The main display shows '0.0' with a 'g' below it.</p>	<p>⇒ Press , the zero display appears. The pictogram →T← is displayed. The pictogram →0← goes out. The weight of the container is now internally saved.</p>
 <p>The display shows a pictogram of a weighing container, a zero with a left arrow, three vertical bars, a full battery icon, and 'CAL'. The main display shows '103' with a 'g' below it.</p>	<p>⇒ Place goods to be weighed in the weighing container. The net weight of the goods to be weighed is displayed.</p>
 <p>The display shows a pictogram of a weighing container, a zero with a left arrow, three vertical bars, a full battery icon, and 'CAL'. The main display shows '- 103' with a 'g' below it.</p>	<p>⇒ The weight of the weighing container will be displayed as a minus number after removing the weighing container.</p>
 <p>The display shows a pictogram of a weighing container, a zero with a left arrow, three vertical bars, a full battery icon, and 'CAL'. The main display shows '0.0' with a 'g' below it.</p>	<p>⇒ The tare weight is saved until it is deleted. Remove the load from the balance and press . The zero display appears, the pictogram →T← goes out and →0← will be displayed again.</p>

8.3 Parts counting

Before the balance can count parts, it must know the average part weight (i.e. reference). Proceed by putting on a certain number of the parts to be counted. The balance determines the total weight and divides it by the number of parts (the so-called reference quantity). Counting is then carried out on the basis of the calculated average piece weight.

i	The higher the reference quantity the higher the counting exactness.
----------	---




Balance display	Operation
	⇒ Start balance by pressing  The balance will carry out a self-test. Wait for „0.0“ display.
	⇒ Press  Appears the reference quantity „10“.
i	By pressing the  key several times, the reference quantities 10, 20, 50, 100, 200 can be set
	⇒ Place as many parts to count on the weighing plate as the set reference quantity requires. ⇒ Confirm by pressing 
 (Example)	The balance is now in parts counting mode and counts all units on the weighing plate.
	⇒ Press  The balance is in weighing mode and displays the weight of the counted parts.
	⇒ Remove the load ⇒ Back to weighing mode by pressing again the  button.

8.4 Adding

It is useful, when a mixture of various components is weighed in a tare container and finally for control purposes the sum weight of all weighed components is required.

Example:

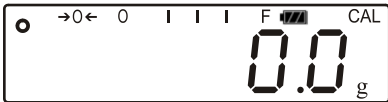





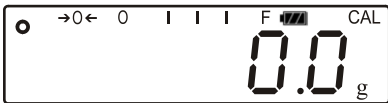

Balance display	Operation
	⇒ Start balance by pressing Whilst the balance is carrying out a selftest, press . „F1 Unt“ appears,
	⇒ Press repeatedly until „F4 ACC“ appears
	⇒ Press , „ACC OFF“ appears
	⇒ Press button to adjust „ACC on“
	⇒ Acknowledge with and press Print to return into weighing mode
	Put on the first sample The weight of the first sample is displayed ⇒ Press
	„ACC 1“ is shortly displayed, then appears once more the weight of the first test item.
	⇒ Remove the first sample ⇒ Put on the second sample The weight of the second sample is displayed. ⇒ Press
	„ACC 2“ is shortly displayed

	<p>Then appears for 2 seconds the weight of both test items, before the weight of the second test item will appear again.</p>
<p>With the further samples proceed as described above.</p>	
<p>In order to display the total weight of the samples, remove weight from the weighing plate, wait for the zero display and press . The number of weighing actions and the total weight of the samples will be displayed shortly.</p>	
<p>i</p>	<p>In order to delete the stored values, press  and while the sum of the test items is displayed, press , the balance will return to „0.0 g“.</p>

8.5 Percent weighings

Percent weighing allows to display weight in percent, in relation to a reference weight.

Percent weighing allows to display weight in percent, in relation to a reference weight.

Balance display	Operation
	⇒ Start balance by pressing  The balance will carry out a self-test. Wait for the „0.0 g“ display.
	⇒ Put a reference weight on the weighing plate, which corresponds to 100%
	⇒ Press  In the display appears „100.00%“
	⇒ Remove reference weight The display returns to „0.0 %“
	⇒ Put on specimen
	In the display appears the percentage of the specimen with reference to the reference weight.
	⇒ Back to weighing mode by pressing again the  button.

9 The menu




9.1 Menu map

In the menu can be made 8 different settings and the adjustment.



9.2 Call up menu



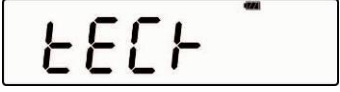




The menu is called-up by pressing the  key, while the balance is carrying out the self test. In the display appears the message „F1 UNT“

9.3 Navigation in the menu

Key	Direction in the menu	Description
	↓	Select menu and pass through menu items from top to bottom
	→	Select current element
	↑	Exit the current element, return to weighing mode

9.4 Menu overview

Balance display	Operation
	Selection of weighing units
	Background lighting on/automatic/off <ul style="list-style-type: none"> • EL on: Background lighting on • EL Au: Background lighting switches on automatically EL off: Background lighting off

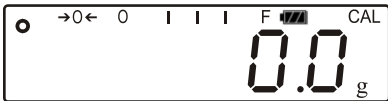







	<p>RS -232 interface S 232:</p> <ul style="list-style-type: none"> • P Prt: connect with printer <ul style="list-style-type: none"> ▪ b XXXX: Baud rate can be selected between 9600, 600, 1200, 2400, 4800 ▪ LP-50: not documented <ul style="list-style-type: none"> • EnG: Selection of Language English • Chi: Selection of Language Chinese ▪ tP: Standard printer setting • P Cont: Continuous printout • P AUto: autom. Printout • P ASk: Data output via remote control commands <p>S USb: not documented</p>	
	<p>Adding (see chap. 8.4): ⇒ ACC of: Adding disabled ⇒ ACC on: Adding enabled</p>	
	<p>Press :</p> <p>Pin: Press ,  and  successively, „P1 Lin“ appears</p>	
	P1 Lin	Linearization (see chapter 6.8)
	P2 CAL	Not documented
	P3 Cont	
	P4 AZn	
	P5 GrA	
	P6 CAP	

9.5 Menu settings


9.5.1 Switch over weighing units (Unit)

The weighing units are switched on or off via the menu.

Activate function:

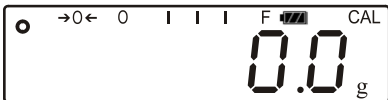









Balance display	Operation
	⇒ Start balance by pressing  As long as the balance carries out a self test, press  .
	„F1 Unt“ is displayed.
 (Example)	⇒ Press  Weighing unit is switched on or off. ⇒ Use  to switch weighing unit on or off ⇒ Using  change to the next unit.

Switch-over weighing unit:

By pressing the  key in the weighing mode it is possible to switch over between the activated units.

9.5.2 Display background illumination

In the menu the display background illumination can be switched on or off.
To achieve this, follow the sequence of operations below:

Balance display	Operation
	⇒ Start balance by pressing  As long as the balance carries out a self test, press  .
	„F1 Unt“ is displayed.
	⇒ Press  „F2 bL“ is displayed.
 (Example)	⇒ Press  „EL on“ or „EL off“ resp. „EL AU“ is displayed ⇒ Press  to switch over between the three settings ⇒ Press  to activate the displayed setting.

Display	Adjustment	Function
„EL on“	Background illumination on	Contrastful display which can also be red in the darkness.
„EL off“	Background illumination off	Battery saving
„EL AU“	The background illumination will be switched off automatically 10 sec after having reached a stable weighing value.	Battery saving

10 Data output RS232

You can print out weighing data via the RS 232C interface.

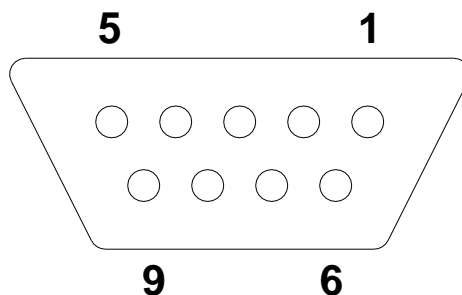
The following conditions must be met to provide successful communication between the weighing balance and the printer.

- Use a suitable cable to connect the weighing balance to the interface of the printer. Faultless operation requires an adequate KERN interface cable.
- Communication parameters (baud rate, bits and parity) of balance and printer must match.

10.1 Technical data

Connection	9 pin d-subminiature bushing
Baud rate	9600
Parity	8 bits, no parity

10.2 Pin allocation of balance output bushing:



Pin 2 input
Pin 3 output
Pin 5 signal earth

10.3 Printer operation - printout examples

Standard printout “weighing data“

GS:	8.65 g	Gross weight
-----	--------	--------------

Printout piece counting

PCS:	10 pcs	Reference quantity
UW:	0.861 g	Individual weight
GS:	8.61 g	Total weight

Printout adding

NO:	1	First weighing
GS:	35.07 g	Weight of first weighing
NO:	2	Second weighing
GS:	8.62 g	Weight of second weighing
NO:	3	Third weighing
GS:	8.00 g	Weight of third weighing
NO:	3	Number weighing processes
GS:	51.69 g	Total of all individual weighings

Printout percent

PERC:	30,19 %	Weight value in percent
-------	---------	-------------------------

10.4 Continuous data output

		,			-/U							k	g	CR	LF
-Header 1-			-Header 2-			-- Weighing data --					Weighing unit		Final character		


Head line 1: ST=stable, US=instable
 Head line 2: NT=net, GS=gross

10.5 Remote control instructions

The remote control commands are sent from the remote control unit to the balance as ASCII code. After the balance having received the s/w/t commands, it will send the following data.

T	Function:	Tare balance
Z	Function:	Set balance to zero
W	Function:	Weighing value for the weight (stable or unstable) is sent via the RS232 interface
S	Function:	Stable weighing value for the weight is sent via the RS232 interface
P	Function:	In counting mode the number of the pieces to be counted is displayed in „Pcs“

11 Error messages

Err 3	Incorrect adjusting weight	Put on correct adjustment weight (see chap. 1; Technical Data)
Err 4	Zero range exceeded	Remove load and press  to reset balance to zero.
Err 5	Keyboard error	Inadmissible input
Err 6	Electronic error	Switch balance off and on again. If the error message remains displayed, please contact your dealer.

12 Service, maintenance, disposal

12.1 Cleaning

Before cleaning, please disconnect the appliance from the operating voltage.

Please do not use aggressive cleaning agents (solvents or similar agents), but a cloth dampened with mild soap suds. Ensure that no liquid penetrates into the device and wipe with a dry soft cloth.

Loose residue sample/powder can be removed carefully with a brush or manual vacuum cleaner.

Spilled weighing goods must be removed immediately.

12.2 Service, maintenance

The appliance may only be opened by trained service technicians who are authorized by KERN.

Before opening, disconnect from power supply.

12.3 Disposal

Disposal of packaging and appliance must be carried out by operator according to valid national or regional law of the location where the appliance is used.

13 Instant help

In case of an error in the program process, briefly turn off the balance and disconnect from power supply. The weighing process must then be restarted from the beginning.

Help:

Fault

Possible cause

The displayed weight does not glow.

- The balance is not switched on.
- The mains supply connection has been interrupted (mains cable not plugged in/faulty).
- Power supply interrupted.
- (Rechargeable) batteries are inserted incorrectly or empty
- No (rechargeable) batteries inserted.

The displayed weight is permanently changing

- Draught/air movement
- Table/floor vibrations
- Weighing plate has contact with other objects.
- Electromagnetic fields / static charging (choose different location/switch off interfering device if possible)

The weighing result is obviously incorrect

- The display of the balance is not at zero
- Adjustment is no longer correct.
- Great fluctuations in temperature.
- Warm-up time was ignored.
- Electromagnetic fields / static charging (choose different location/switch off interfering device if possible)

Should other error messages occur, switch balance off and then on again. If the error message remains inform manufacturer.